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Dear Customer,

Thank you for choosing a high-quality LaCont environmental technology product.

To ensure that the entire system functions perfectly, please read the following operating instructions carefully.

This applies to the following models:

- **CHS – PSM – UWS with hinged doors**
- **CHS with sliding doors**

We cannot accept any liability if these instructions are not followed.

We expressly reserve the right to make technical changes that serve to improve our product or that are necessitated by changes in the law, even without prior notice.

These operating/installation instructions may only be reprinted or reproduced, even in part, with the written permission of Lacont Umwelttechnik GmbH.

LaCont Umwelttechnik GmbH
39435 Egel

Tel.: +49 (0)39 268 / 98 96 - 0, Fax: +49 (0)39 268 / 98 96 - 29,

Email: info@lacont.de , Website: www.lacont.de

Table of contents – Operating instructions

1. General safety instructions / Warranty
2. Intended use
3. Technical data / Drawings
4. Transport
5. Installation
6. Commissioning
7. Interior fittings
8. Storage quantities / Combined storage
9. Ventilation
10. Labeling
11. Maintenance / Servicing
12. Disposal

1. General safety information / Warranty

- The cabinets must be set up, operated, and maintained in such a way that the safety of employees is guaranteed against harmful vapors.

- The storage of flammable liquids, aggressive substances, or substances that could cause fire or explosions due to their self-igniting or unstable properties is not permitted.

- When operating your chemical cabinet, please observe the applicable national rules, regulations, and laws that apply in this regard.

- Before commissioning, a functional test must be carried out to detect any damage caused by transport, handling, and improper installation!

- The warranty/guarantee is provided in accordance with our General Terms and Conditions. See also point 12.

- Please observe your duty of care—damage must be repaired immediately and leaks must be removed!

- The door opening area must be kept clear and clean.

- Doors and drawers should only be kept open for loading and unloading – keep doors/drawers closed!

- Please observe the maximum storage quantities, load capacity of the shelves, etc. in accordance with Appendix 1 or the labeling on the door/drawer.

- The chemical cabinets are not intended for storing aggressive substances (e.g., acids and alkalis, corrosive gases) – otherwise there is a risk of corrosion of the interior fittings and the exhaust air components! Please use our special acid and alkali cabinets for this purpose.

- Before storing the goods (e.g., chemicals), check the resistance of the cabinet surface and the resistance of the drip tray.

- The collection device (floor tray) must be able to collect 10% of the total quantity stored above it, but at least the contents of the largest single container. In water protection areas, 100% of the stored quantity must be collected, provided that storage is permitted. (valid for Germany)

- In addition, the applicable national regulations must be observed and complied with.

- When handling hazardous substances or toxins, suitable personal protective equipment must be worn!

2. Intended use

storeLAB® chemical cabinets are used for the storage of water-polluting and/or toxic substances in sealed containers in workrooms, provided that storage is compatible with the protection of employees.

3. Technical data / drawings

The technical data and drawings for the individual model variants can be found in Appendix 1 starting on page 40.

3.1 Construction

Body: Powder-coated sheet steel

Features: Height-adjustable insert trays or tray bases, shelves, and pull-out trays

Collection tray in accordance with steel tray guidelines with Ü mark (not for CHS Eco)

4. Transport

The storeLAB® chemical cabinet may only be lifted from the bottom using suitable means of transport and must be transported in an upright position. It is essential to avoid tilting during lifting and transport. Always lift the cabinet from the side or rear. To avoid damage to the doors, never lift it from the door side. The lifting surface of the hand truck, pallet truck, or similar must be at least the same size as the base of the safety cabinet.

CAUTION!

Risk of crushing – use of PPE

5. Installation

The floor must be level in order to support the weight of the storeLAB® chemical cabinets when fully loaded.

To prevent tipping and ensure proper use of the doors, storeLAB® chemical cabinets must be installed horizontally.

This is the only way to ensure that the cabinets cannot tip over.

Due to the location of the actual place of use/operation, it must be ensured that the chemical cabinets

- cannot be damaged by vehicles.
- not be exposed to any direct or indirect heat sources - be protected from moisture

When handling harmful or toxic substances, wear suitable personal protective equipment!

The location of the exhaust air connections can be found in Appendix 1 (see sketches) starting on page 40.

6. Commissioning

- 6.1 The storeLAB® chemical cabinet must be installed, operated, and maintained in such a way that the protection of employees is guaranteed.
- 6.2 The storeLAB® environmental wall cabinet must be installed, operated, and maintained in such a way that the safety of employees is guaranteed.
 - The installation instructions can be found in the appendix on page 15.

ATTENTION!

When handling harmful or toxic substances, suitable personal protective equipment must be worn!

7. Interior

7.1 Insert tray / shelves

The insert trays and shelves (CHS with sliding doors) supplied are factory-fitted in the cabinet. To position the insert tray, lift the tray and rotate it around its horizontal axis until it can be pulled out of the cabinet. Then lift the shelf support and remove it from the mounting bracket. Assembly is carried out in reverse order.

CAUTION! Risk of crushing if handled improperly - use PPE

7.2 Pull-out trays

The pull-out trays supplied and the pull-outs with tray supports are installed in the cabinet at the factory. To position the pull-out tray, pull it forward, release the lock on the pull-out and remove it horizontally from the cabinet. Pull the pull-out supports with pull-out from the holes in the adjustment rails and turn them 90° upwards. Pull the pull-out support, which is now horizontal, out of the holes in the adjustment rails. Installation in the new position is carried out in the reverse order of the procedure described above.

8. Storage quantities / Combined storage

The applicable national regulations must be observed.

For maximum storage quantities, see Appendix 1 from page 44 onwards!

9. Ventilation

PSM / UWS / CHS Eco cabinets do not have an exhaust air connection; all other chemical cabinets have an exhaust air connection (Ø80 mm). Please observe the applicable national rules, regulations, and laws regarding the prescribed air exchange and exhaust air ducting. The location of the exhaust air connections can be found in Appendix 1. The storeLAB® chemical cabinets can also be operated without technical exhaust air.

The applicable national rules, regulations, and laws must be observed in this regard.

10. Labeling

Labeling according to TRGS / ISO 3864 / GHS-VO.

11. Maintenance

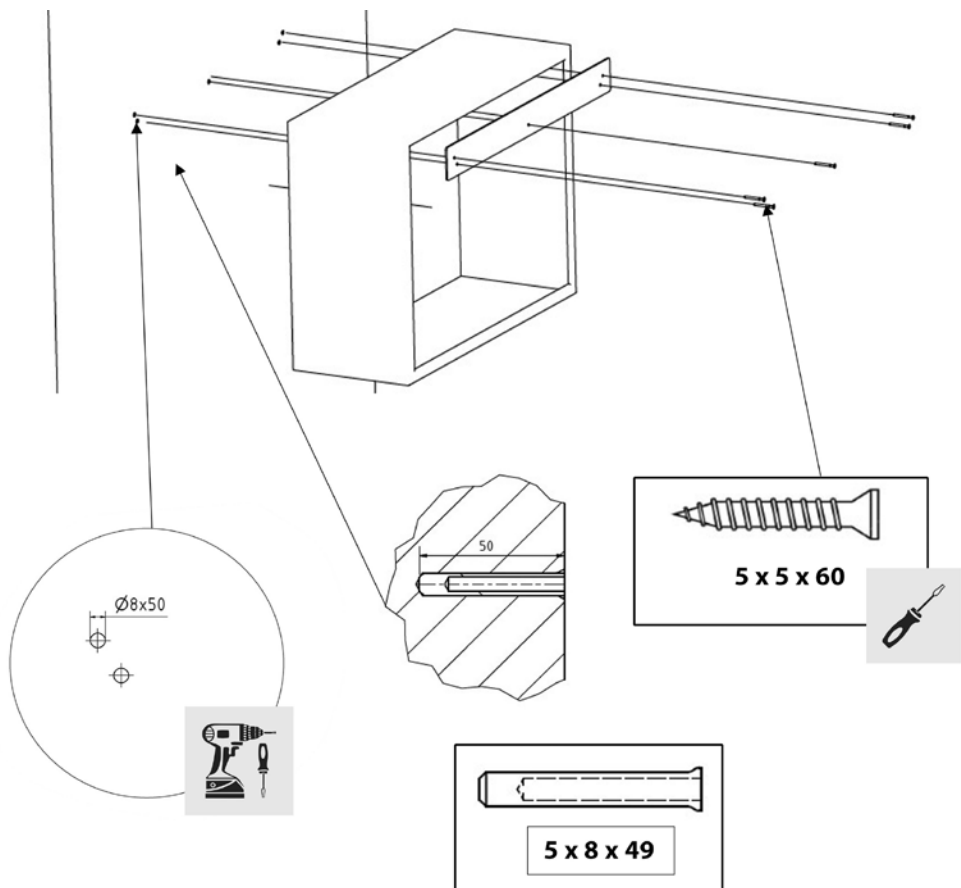
11.1 Daily inspection

Spilled liquids must be cleaned up immediately and disposed of properly. Check the cabinet regularly for mechanical damage, contamination, and corrosion.

12. Disposal

The storeLab® chemical cabinet can be completely dismantled. The individual components can be recycled separately. The applicable national and local disposal regulations must be observed. To protect resources, parts of the chemical cabinet or the entire cabinet should not be disposed of in bulky waste or household waste.

Assembly instructions for installation Environment Wall cabinets

5 x 8 x 49 dowels and 5 x 5 x 60 screws are included. Check the condition of the wall.

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storeLAB® Chemicals-Cabinet (*included GL-Version*)
- Cabinets for Pesticides with hinged door
- Environmental cabinet with hinged door
- Chemical cabinet with sliding door

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to which this declaration relates, is in conformity with the relevant provisions of the following standards or other normative documents.

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**EG-Maschinenrichtlinie 2006/42/EG, L'157/65
vom 9.6.2006**
EEC-Machinery Directive 2006/42/EEC

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Applied national
standards (germany)

DIN 4844-1 Safety colours and safety signs

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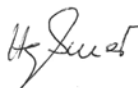
The operation company is responsible for installing the products in accordance to national regulations, which may apply in respect to the installation, use and also for regular inspections.

Egeln, den

01.07.2025

Datum

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D - 39435 Egelu / GERMANY

Phone: +49 (0) 39 268 / 98 96 - 0

Fax: +49 (0) 39 268 / 98 96 - 29

E-Mail: info@lacont.de

Internet: www.lacont.de